


MWS-Precisa-Application Laboratory

Auerstrasse 17
CH 9435 Heerbrugg
e-mail: monika.blum@precisa.ch

Tel.: ++ 41 71 727 13 80
Fax: ++41 71 727 13 99

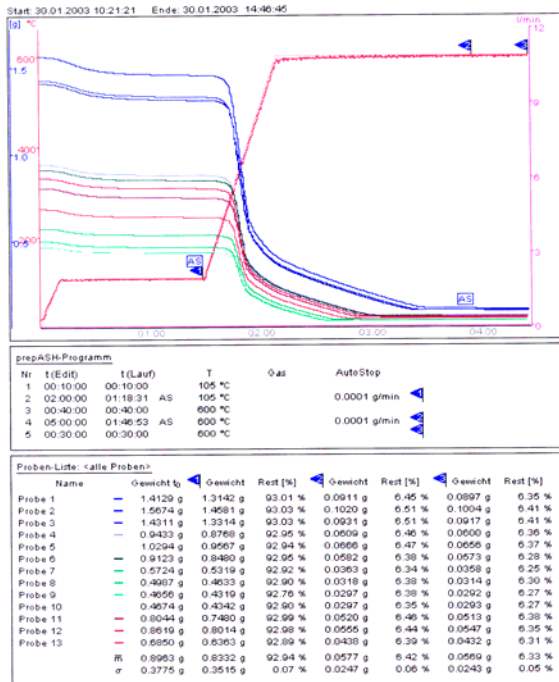

ASH in PP/GF recycled material

All fundamental data and operating parameters, such as changes in sample weight, temperature, gases, etc. are stored and graphically displayed (see graphic example on the left)
Automatic calculation of results.

Methode:

Step	time	temperature 1	temperature 2	Auto Stop	gas
1	45 min	25 °C	900 °C		
2	3 h	900C	900°C		

Result:

Sample amount	2 g
Ash (Glass fiber content)	27.18 %
Standard deviation	0.19 %
n	29

Heerbrugg, 22.10.2003

Monika Blum